A BETTER VIEW OF THE WORLD

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EUROPEAN
OPTICS
since
1933

MEASUREMENT OF OPTICAL PARAMETERS OF MICROSCOPE OBJECTIVES

Content

- **WHAT IS A MICROSCOPE OBJECTIVE (MO)**
- **WHAT WE CAN KNOW ABOUT MO**
- **WHAT CAN BE MEASURED**
- **WHAT CAN BE CUSTOMISED**

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
MEASUREMENT OF OPTICAL PARAMETERS OF MICROSCOPE OBJECTIVES

What is a microscope objective?

At its simplest definition: mo is a very high-powered **magnifying lens**, with very short **focal length**.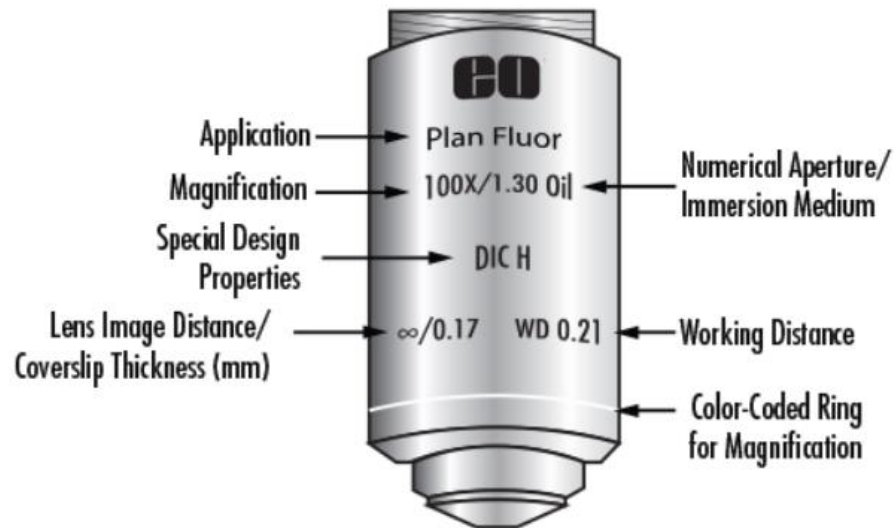


MEASUREMENT OF OPTICAL PARAMETERS OF MICROSCOPE OBJECTIVES

How to sort the Microscope objectives?

MICROSCOPE OBJECTIVES CAN BE SORTED BY:

- Magnification (1x, 10x, 20x, 100x, ...)
- Numerical aperture (resolution) (0,2, 0,4, 0,7, 0,9, 1,2, ...)
- Manufacturer (nikon, zeiss, olympus, leica, mitutoyo, ...)
- Working distance (short WD, long WD, ...)
- Immersion or cover glass correction (yes / no)
- Application (aberration correction type)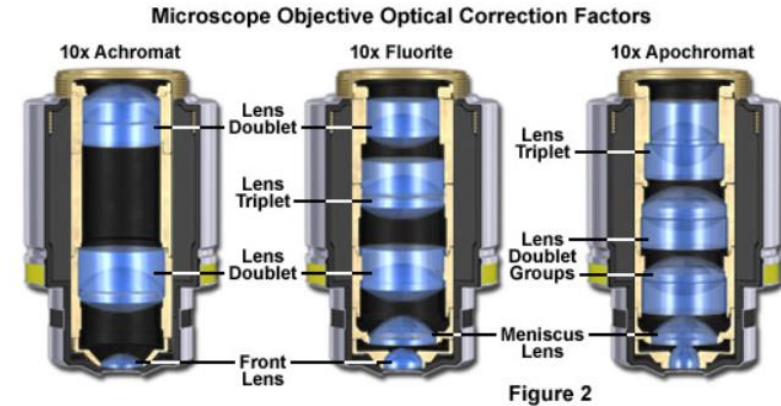

Microscope Objective Correction for Optical Aberration

| Objective Specification | Spherical Aberration | Chromatic Aberration | Field Curvature |
|-------------------------|----------------------|----------------------|-----------------|
| Achromat | 1 Color | 2 Colors | No |
| Plan Achromat | 1 Color | 2 Colors | Yes |
| Fluorite | 2-3 Colors | 2-3 Colors | No |
| Plan Fluorite | 3-4 Colors | 2-4 Colors | Yes |
| Plan Apochromat | 3-4 Colors | 4-5 Colors | Yes |


MEASUREMENT OF OPTICAL PARAMETERS OF MICROSCOPE OBJECTIVES

What are the parameters published by supplier?

SPECIFICATION – Typical published objectives parameters

| Specifications Learn More Accessories Technical Information Related Products | | | |
|--|------------------------|--|--|
| Magnification: | 40X | Numerical Aperture NA: | 0.6 |
| Compatible Tube Lens Focal Length (mm): | Focal Length: 164.50mm | Compatible Cover Glass Thickness (mm): | N/A |
| Depth of Field (µm): | 0.76 | Field Number (mm): | 23 |
| Focal Length FL (mm): | 4.10 | Length excluding Threads (mm): | 42.8 |
| Manufacturer: | ZEISS | Maximum Diameter (mm): | 30 |
| Model Number: | 422060-9901-000 | Mounting Threads: | M27 x 0.75 |
| Parfocal Length (mm): | 45 | Resolving Power (µm): | 0.46 |
| Style: | Infinity Corrected | Type: | Microscope Objective |
| Weight (g): | 180 | Working Distance (mm): | 2.2 |
| Wavelength Range (nm): | 480 - 650 | Immersion Liquid: | N/A |
| Operating Temperature (°C): | +10 to +40 | Typical Applications: | Fluorescence, BrightField, Differential Interference Contrast, Reflected Light DIC, DIC with circular polarized light, total interference contrast |

Built-in (instalation) parameters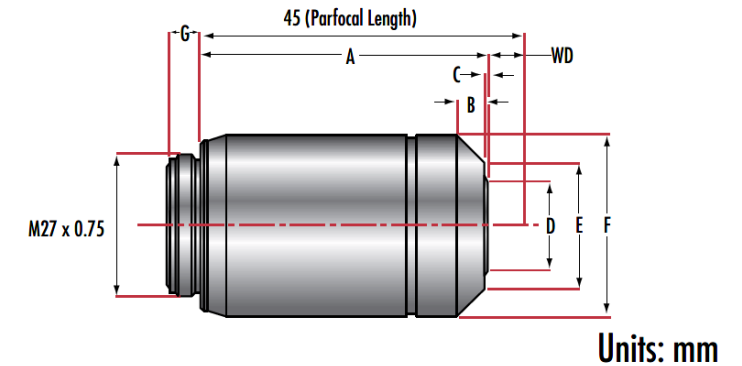




Technical Information

| Stock No. | EC Epiplan | Dimensions (mm) | | | | | | | | W.D. | PFL | Mounting Threads |
|-----------|------------|-----------------|------|------|------|------|------|-----|-------|-------|----------|------------------|
| | | A | B | C | D | E | F | G | | | | |
| #13-822 | 5X | 33.20 | 2.10 | 6.00 | 11.6 | 22.5 | 30.0 | 3.9 | 11.80 | 45.00 | M27x0.75 | |
| #13-823 | 10X | 34.00 | 2.10 | 6.80 | 11.6 | 22.5 | 30.0 | 3.9 | 11.00 | 45.00 | M27x0.75 | |
| #13-824 | 20X | 41.80 | 7.40 | 3.80 | 12.6 | 22.5 | 30.0 | 3.9 | 3.20 | 45.00 | M27x0.75 | |
| #13-825 | 40X | 42.80 | 9.40 | 2.80 | 9.1 | 14.5 | 30.0 | 3.9 | 2.20 | 45.00 | M27x0.75 | |
| #13-826 | 50X | 44.05 | 8.10 | 5.35 | 9.0 | 22.5 | 30.0 | 3.9 | 1.00 | 45.00 | M27x0.75 | |
| #13-827 | 100X | 44.15 | 8.10 | 5.45 | 9.0 | 22.5 | 30.0 | 3.9 | 0.87 | 45.00 | M27x0.75 | |

MEASUREMENT OF OPTICAL PARAMETERS OF MICROSCOPE OBJECTIVES

What about quality?

There does not exist any information about quality in any of manufacturer (MTF, WFE, resolution)? Why?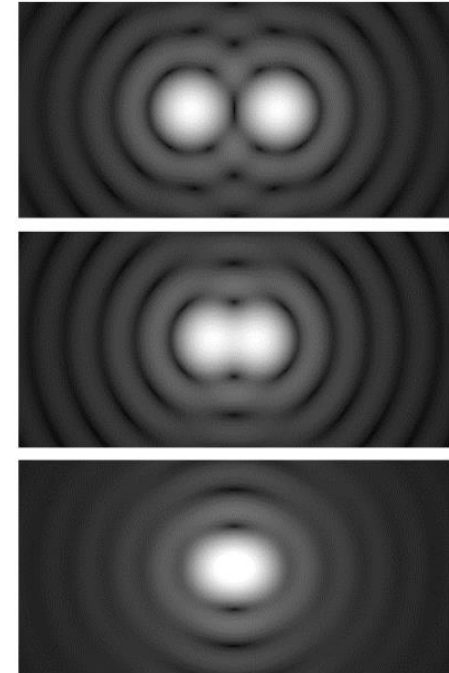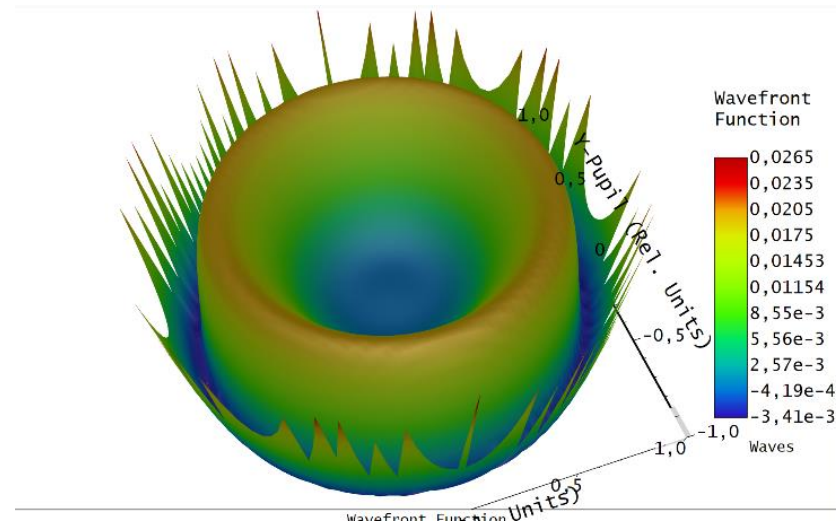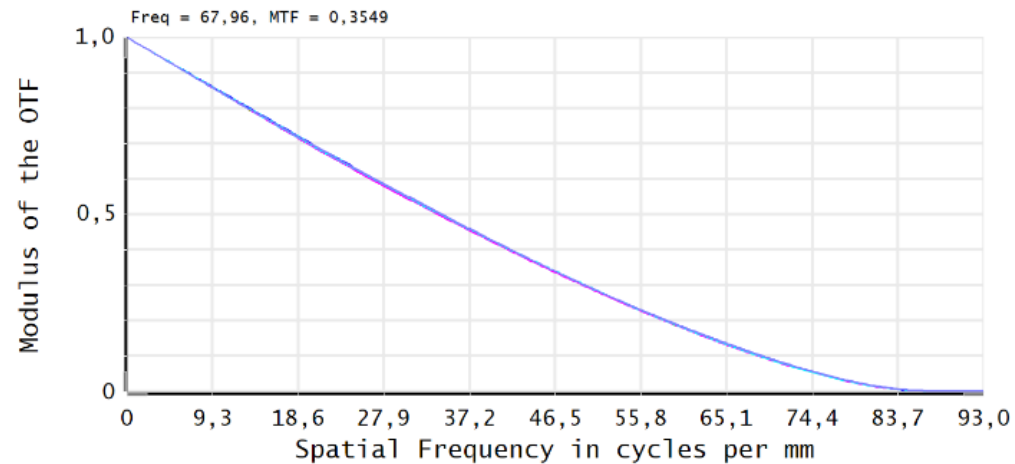


MEASUREMENT OF OPTICAL PARAMETERS OF MICROSCOPE OBJECTIVES

What would we really like to know about the microscope objective?



GEOMETRICAL OPTICAL PARAMETERS:

- Distortion
- Field curvature
- Lateral & longitudinal chromatic aberration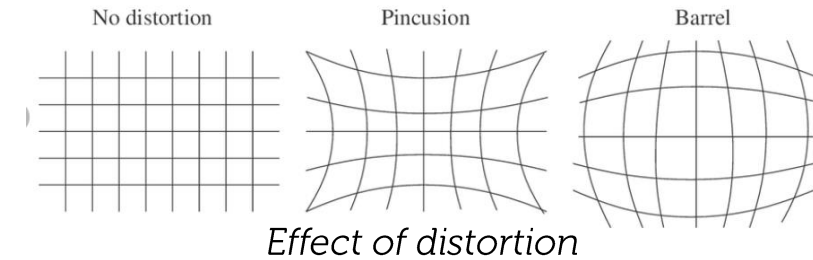

QUALITATIVE OPTICAL PARAMETERS:

- WFE / MTF / resolution

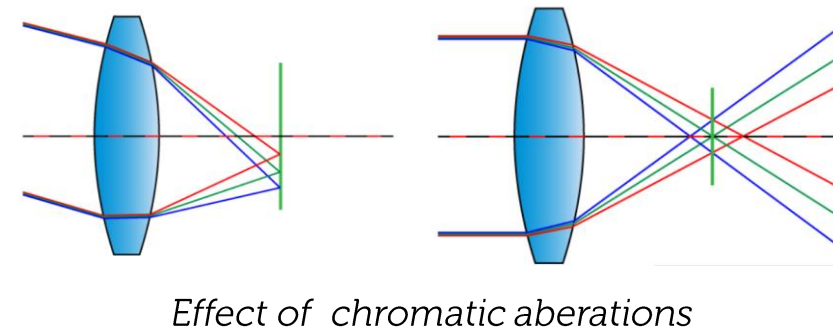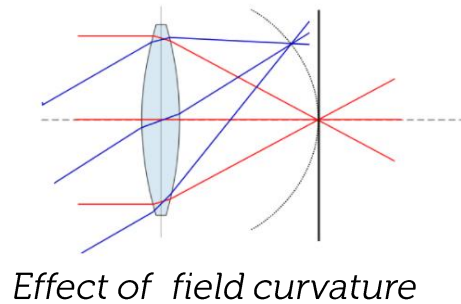MEASUREMENT OF OPTICAL PARAMETERS OF MICROSCOPE OBJECTIVES

A BETTER
VIEW
OF THE
WORLD

How to determine microscope objective quality?

WHAT OPTIONS WE HAVE?

- Ask supplier *
- Hope for good results / condition
- **Measure**

* The only information which is (sometimes) possible to ask from supplier can be a blackbox design (zemax, codev)


- > BUT even if so, this shows just the nominal design
 - > no real system behaviour (after tolerancing)
 - > no volume statistics

MEASUREMENT OF OPTICAL PARAMETERS OF MICROSCOPE OBJECTIVES


A BETTER
VIEW
OF THE
WORLD

How to measure microscope objectives?

Meopta R&D labs has dedicated experimental setups to measure geometrical & qualitative parameters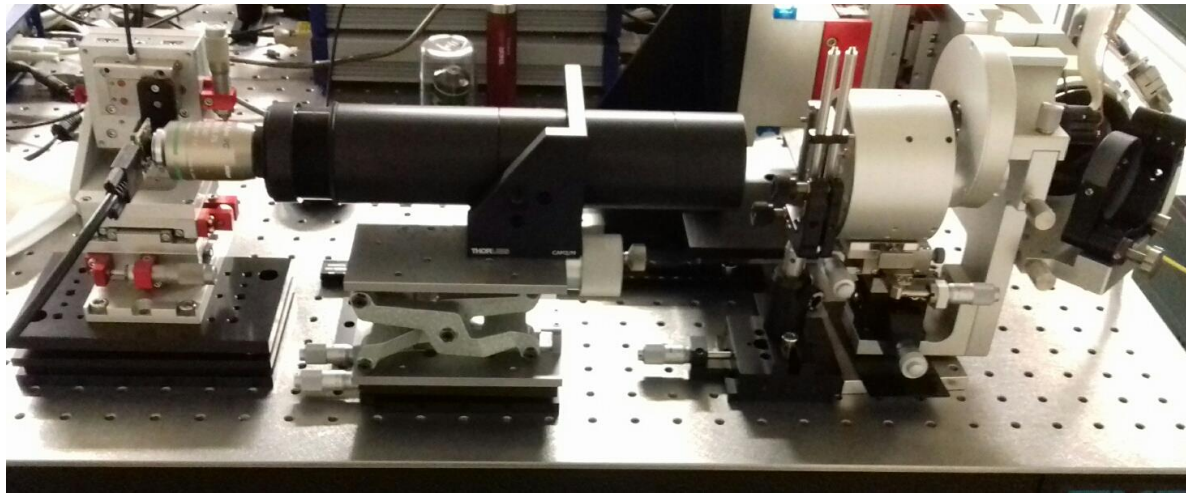

Experimental setup for geometrical parameters measurement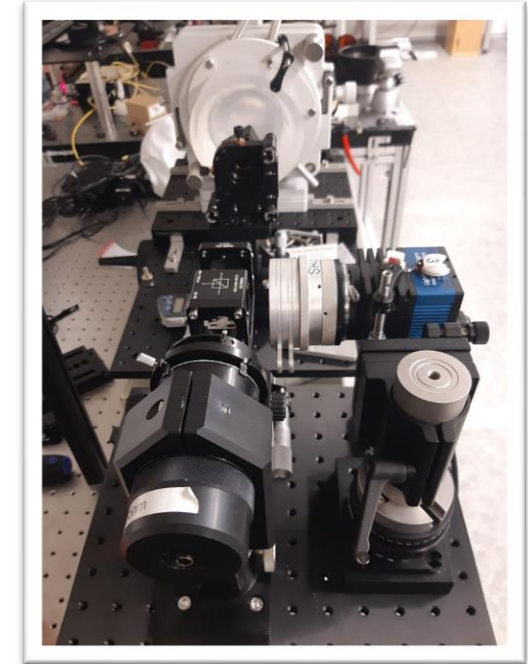




Experimental setup for qualitative parameters measurement

MEASUREMENT OF OPTICAL PARAMETERS OF MICROSCOPE OBJECTIVES

Example of MO measurement / selection

Target: to select the optimal option of MO for specific design / application

| MO selection parameters | |
|-------------------------|-------------|
| NA | 0,7-0,8 |
| Magnification | 20x |
| Spectrum | 350 – 650nm |
| WD | 0,5 – 1,5mm |

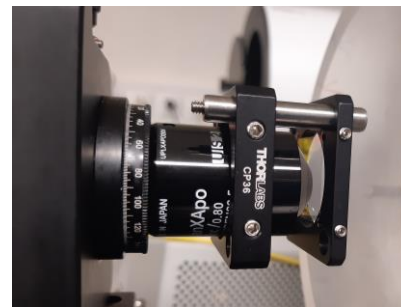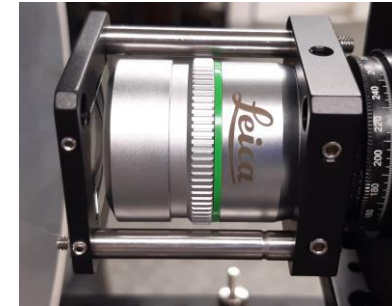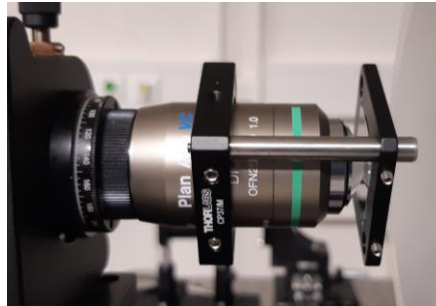
MEASUREMENT OF OPTICAL PARAMETERS OF MICROSCOPE OBJECTIVES

How to measure microscope objectives?

GEOMETRICAL PARAMETERS

Distortion & field curvature results – displacement in lateral & longitudinal axis

| Distortion | |
|------------|-----|
| PV [nm] | |
| Leica | 901 |
| Zeiss | 326 |
| Nikon | 517 |
| Olympus | 329 |


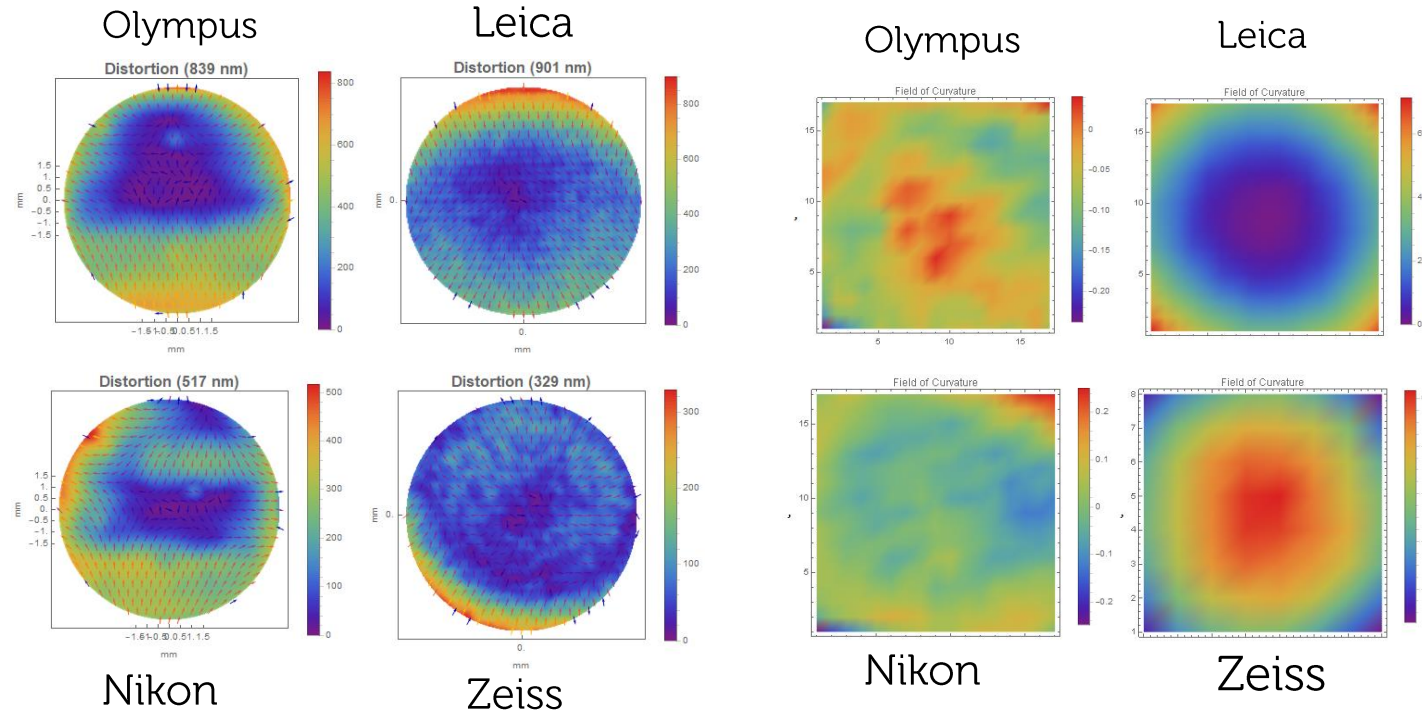
| Field curvature | |
|-----------------|-----|
| PV [nm] | |
| Leica | 750 |
| Zeiss | 430 |
| Nikon | 80 |
| Olympus | 280 |

MEASUREMENT OF OPTICAL PARAMETERS OF MICROSCOPE OBJECTIVES

How to measure microscope objectives?

GEOMETRICAL PARAMETERS



Longitudinal and transversal chromatic aberrations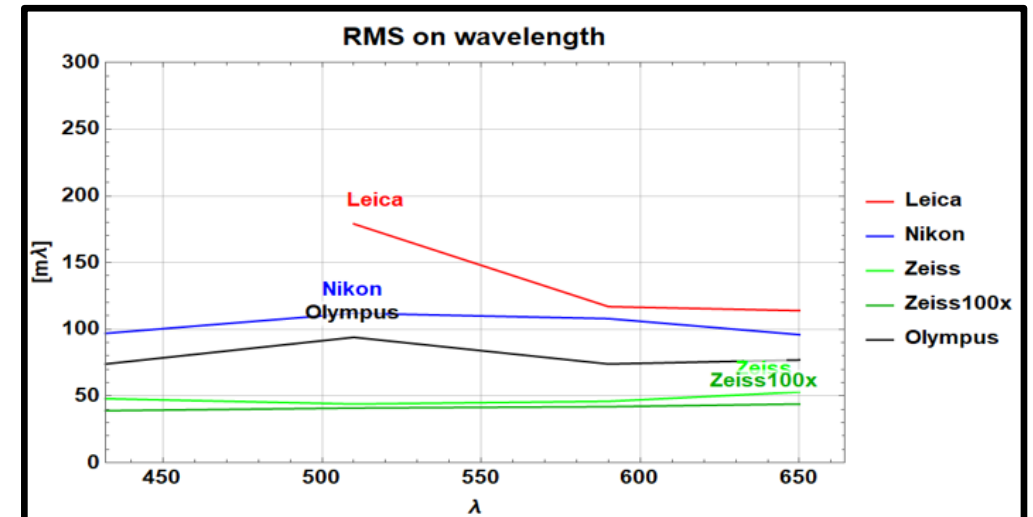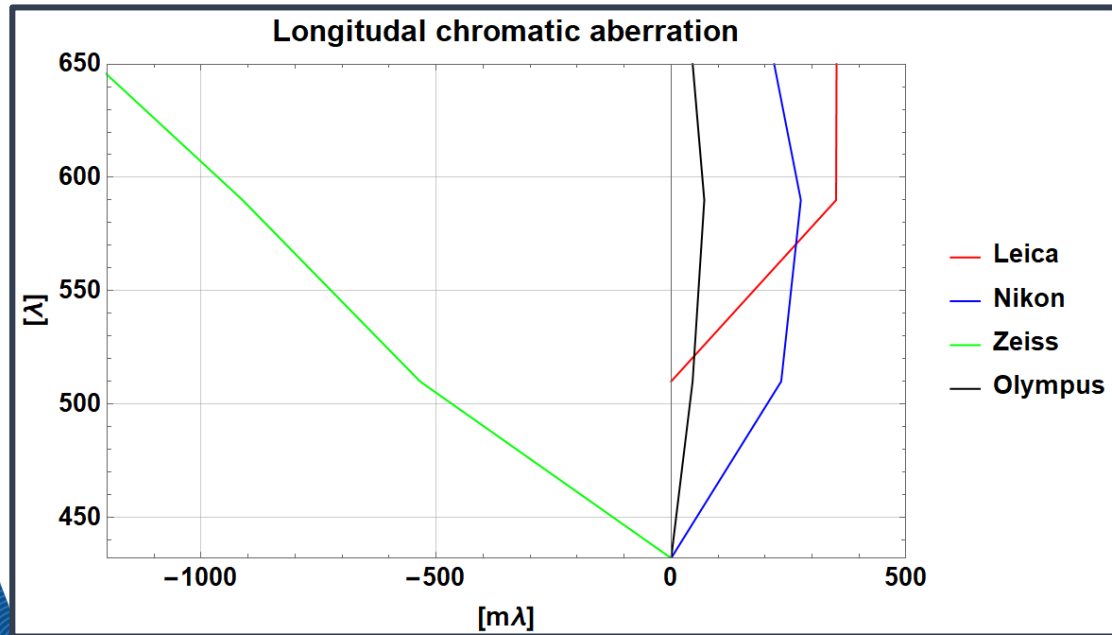

MEASUREMENT OF OPTICAL PARAMETERS OF MICROSCOPE OBJECTIVES

How to measure microscope objectives?

QUALITATIVE PARAMETERS - WFE

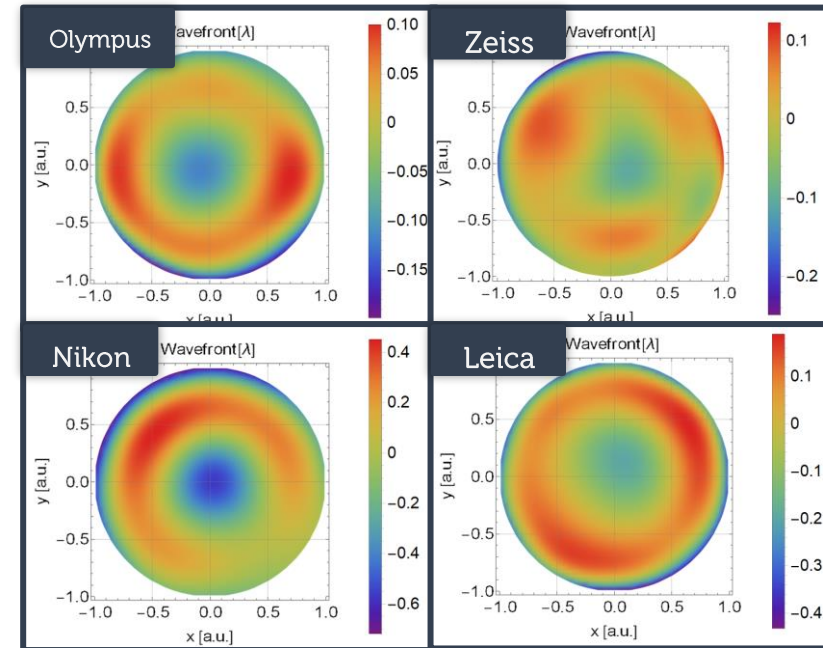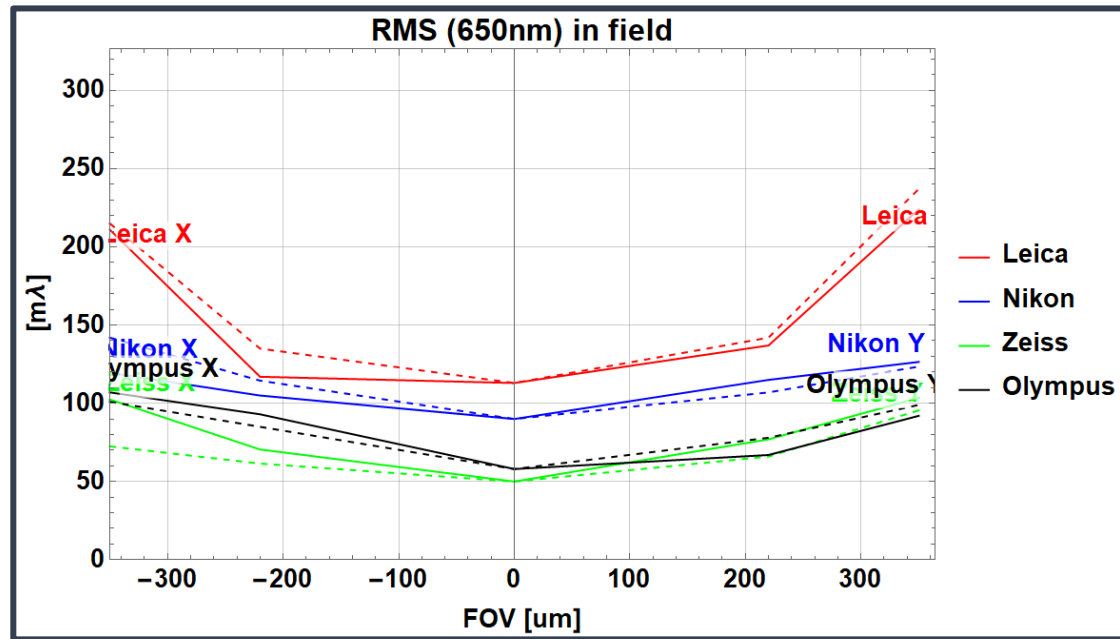| Objektiv | průměr 440um [mλ] | průměr 700um [mλ] |
|----------|-------------------|-------------------|
| Leica | 132.75 | 221.75 |
| Nikon | 110.37 | 127.88 |
| Zeiss | 68.75 | 93.5 |
| Olympus | 80.75 | 99.75 |

MEASUREMENT OF OPTICAL PARAMETERS OF MICROSCOPE OBJECTIVES

How to measure microscope objectives?

Final selection done as a result of technical (quality & geometrical) parameters as well as non-technical parameters (application, space, etc.)

| Result | | | | |
|----------------|---------------|-----------------|---------------|----------------|
| | Score - tech. | Score - non-tec | Complex score | Complex rating |
| Zeiss | 15 | 6 | 21 | 1 |
| Olympus | 11 | 3 | 14 | 3 |
| Nikon | 10 | 5 | 15 | 2 |
| Leica | 1 | 1 | 2 | 4 |

| Technical parameters | | | Non-technical parameters | | |
|----------------------------|--------|--------|--|--------------|--------|
| Wavefront | | | Focal length / Working distance | | |
| | Rating | Points | | Rating | Points |
| Zeiss | 1 | 9 | Zeiss | 8,25 (WD1,3) | 2 |
| Olympus | 2 | 6 | Olympus | 9 (WD0,6) | 0 |
| Nikon | 3 | 3 | Nikon | 10 (WD1) | 1 |
| Leica | 4 | 0 | Leica | 10 (WD0,6) | 0 |
| Geometrical | | | Cover glass | | |
| Distorsion | | | | Rating | Points |
| Zeiss | 1 | 3 | Zeiss | No | 2 |
| Olympus | 3 | 1 | Olympus | Yes | 0 |
| Nikon | 2 | 2 | Nikon | Yes | 0 |
| Leica | 4 | 0 | Leica | Yes | 0 |
| Field curvature | | | Vacuum compatibility | | |
| | | | | Rating | Points |
| Zeiss | 2 | 2 | Zeiss | Yes | 2 |
| Olympus | 3 | 1 | Olympus | Maybe | 1 |
| Nikon | 1 | 3 | Nikon | Maybe | 1 |
| Leica | 4 | 0 | Leica | No | 0 |
| Chromatic aberation | | | Price | | |
| | Nedef | | | Rating | Points |
| Zeiss | | 1 | Zeiss | | 0 |
| Olympus | 1 | 3 | Olympus | | 2 |
| Nikon | 2 | 2 | Nikon | | 3 |
| Leica | 3 | 1 | Leica | | 1 |

MEASUREMENT OF OPTICAL PARAMETERS OF MICROSCOPE OBJECTIVES

A BETTER
VIEW
OF THE
WORLD

What is the other option?

Meopta designs a customised microscope objectives and microscope systems for customers from semiconductor, FPD, fluorescence microscopy, medical, etc.




MEASUREMENT OF OPTICAL PARAMETERS OF MICROSCOPE OBJECTIVES

A BETTER
VIEW
OF THE
WORLD

Example of meopta design?

| | nominal design | units |
|------------------------|----------------|--------|
| Wavelength range | 435 - 485 | |
| Object NA | 0,4 | - |
| Image NA | 0,022 | - |
| Object FOV (semi-diam) | 4 | mm |
| Image FOV (semi-diam) | 80 | mm |
| Paraxial magnification | 20 | - |
| WFE | 0,008 | mλ RMS |
| Strehl ratio | 0,997 | - |

Example of customized microscope systems specs

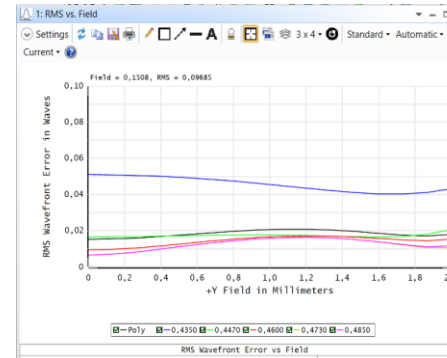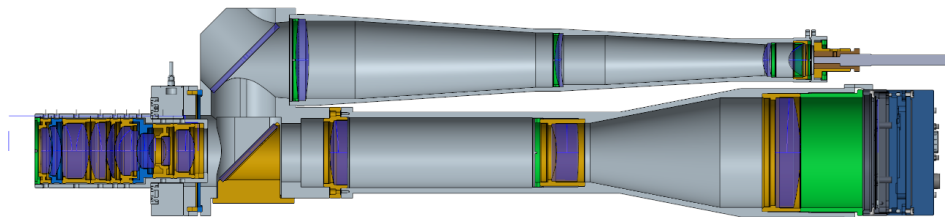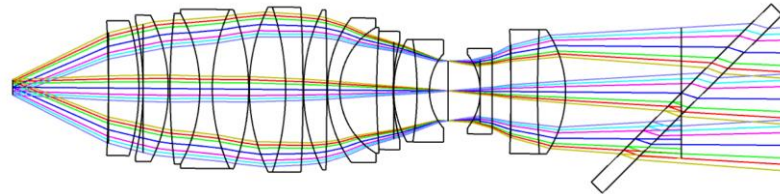MEASUREMENT OF OPTICAL PARAMETERS OF MICROSCOPE OBJECTIVES

Summary

- MICROSCOPE OBJECTIVES (MO) GENERALLY
- WHAT THE MANUFACTURERS PUBLISH ABOUT MICROSCOPE OBJECTIVES (WHY THERE IS NO INFO ABOUT QUALITY)
- WHAT CAN BE MEASURED + EXAMPLE OF MO SELECTION
- EXAMPLE OF CUSTOMISED MO

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